

LD

AUTOCLAVES VERTICALES



Série **LD** d'autoclaves verticales de grand volume pour stérilisation à vapeur de pression en acier inoxydable

Modèle : **LD -100** \ **LD-150** \ **LD-200**

Volume : 100L \ 150L \ 200L

Matière: acier inoxydable complet

- ◇ Structure sur roues pour un facile déplacement et ouverture du couvercle rapide
- ◇ contrôle automatique et programme cyclique antiseptique
- ◇ écran LCD indicateur de l'état de travail
- ◇ verrouillage de sécurité en présence de pression
- ◇ champ de température 50 °C -134 °C
- ◇ sélection des temps de stérilisation de temps :0-99h
- ◇ valve de sécurité à déclenchement automatique pour surpression 0,145-0.208Mpa
- ◇ Arrêt automatique avec signal sonore à la fin du cycle de stérilisation
- ◇ Avec protection contre manque ou insuffisance d'eau
- ◇ Complètement réalisée en acier inoxydable
- ◇ Manomètre de contrôle de la pression/température



Tableau de référence pour les articles destinés à être stérilisés

Articles à stériliser	Temps de stérilisation à la température programmée (min)	Température de stérilisation (°C)	Pression relative (MPa)
Caoutchouc	15	~121	0.1~0.105
Tissus	30-45	121~126	0.105~0.14
Verrerie	15	121~126	0.105~0.14
Instruments	10	121~126	0.105~0.14
Solutions	20-40	121~126	0.105~0.14
Vaisselle	20	121~126	0.105~0.14

Model	Volume	Alimentation	Consommation	Dimensions de la chambre	Poids	Dimensions
LD-100	100L	380V	6,0 Kw	Φ 500×550 mm	100kg	700×760×1320 mm
LD-150	150L	380V	7,5 Kw	Φ 500×760 mm	120kg	730×760×1480 mm
LD-200	200L	380V	9.0 Kw	Φ 500×1000 mm	140kg	730×780×1690mm



ESSE 3 Via Garibaldi 30
14022Castelnuovo D.B. (AT)
tel +39 011 99 27 706
fax +39 011 99 27 506
e-mail esse3@chierinet.it
web : www.esse3-medical.com



LD VERTICAL AUTOCLAVES

Series Vertical Type Large-volume Stainless Steel Pressure Steam Sterilizer

Model : **LD -100** \ **LD -150** \ **LD -200**

Volume : **100L** \ **150L** \ **200L**

Material : Full stainless steel

- ◇ Hand wheel type of quick-open door structure
- ◇ Automatic control antiseptic cyclic program
- ◇ LCD screen indicate working state
- ◇ Pressure safety lock
- ◇ Temperature scope (50°C-134°C)
- ◇ Time scope (0-99h)
- ◇ Overpressure auto-discharging 0.145-0.208Mpa
- ◇ Automatic shut down with beep reminding after sterilization
- ◇ With water break protecting equipment
- ◇ Full stainless steel

Reference Table for Articles to Be Sterilized

Articles to Be Sterilized	Temperature Retaining Time (min)	Relative Temperature for Steam (°C)	Relative Pressure for Steam (MPa)
Rubber	15	~121	0.1~0.105
Dressings	30-45	121~126	0.105~0.14
Glassware	15	121~126	0.105~0.14
Devices & Instruments	10	121~126	0.105~0.14
Solutions	20-40	121~126	0.105~0.14
Tableware	20	121~126	0.105~0.14

Model	Volume	Power Source	Consumed Power	Dimension of sterilizing chamber	Weight	Dimension
LD-100	100L	380V	6,0 Kw	Φ500×550 mm	100kg	700×760×1320 mm
LD-150	150L	380V	7,5 Kw	Φ500×760 mm	120kg	730×760×1480 mm
LD-200	200L	380V	9.0 Kw	Φ500×1000 mm	140kg	730×780×1690mm



ESSE 3 Via Garibaldi 30
14022Castelnuovo D.B. (AT)
tel +39 011 99 27 706
fax +39 011 99 27 506
e-mail esse3@chierinet.it
web : www.esse3-medical.com

